Moticam A Series

Quick Start Guide

Start your discovery journey with your new Moticam

ALL IN ONE BOX

Accessories included in the package:

1. Moticam

2. C-ring

3. Focusable lens (12mm)

4. Eyepiece adapters (2)

5. USB cable

6. Macro-tube

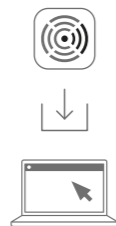
7. Calibration Slide

8. Software (download link)


<https://ws.motic.com/mo1/moticonnect>

1

Download the MotiConnect Software from:
<https://ws.motic.com/mo1/moticonnect>
and install it in your computer.

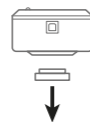2


Remove the Dust Cap from the camera.
To find out how to use your Moticam, follow the steps below according to your needs:
3A - On a trinocular microscope
3B - Attached to the eyepiece
3C - Without a microscope as a Macro camera

Continue to steps 3A, 3B or 3C


3A

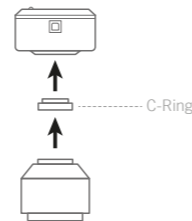Using the camera on a **Trinocular Microscope**

To use your Moticam on a **trinocular microscope**, please attach the C-Mount adapter to the camera.
**Please note that, depending on the type of adapter and microscope, the C-Ring may be necessary to properly connect the camera and C-Mount adapter.*

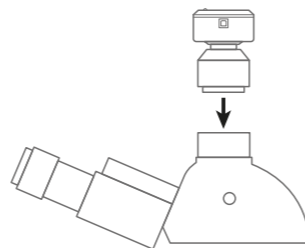4A


Attach the combination of camera and adapter to the trinocular tube of your microscope.

Continue to step 5


3B

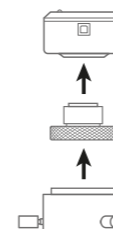Using the camera **on an eyepiece**

To use your Moticam with an **eyepiece adapter**, please attach the focusable lens to the camera. Choose the right size eyepiece adapter, then proceed to connect it to the focusable lens.

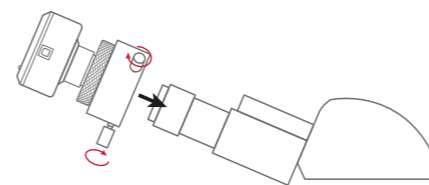4B


Attach this assembly to your microscope's eyepiece by tightening the centering screws on the eyepiece adapter.

Continue to step 5

3C

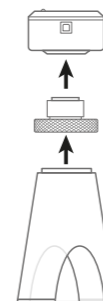Using the camera as a **Macro Camera**

To use the Moticam as a **Macro Camera**, please attach the focusable lens to the camera, then proceed to connect the Macro Tube onto the lens.

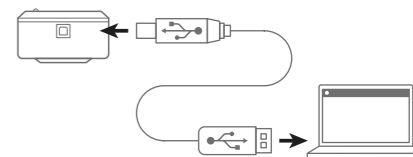
4C

Place this assembly over the item you wish to investigate and ensure there is adequate lighting. (If necessary, a desk lamp is a good solution).

Continue to step 5

5

Connect the USB cable to the camera and then plug it into your computer.

6

Open the MotiConnect Software, click on "Camera Device" to select your camera model, and see your live image.

Don't miss our video-tutorials:

MotiConnect
FIRST STEPS

MotiConnect
CALIBRATION &
MEASUREMENTS

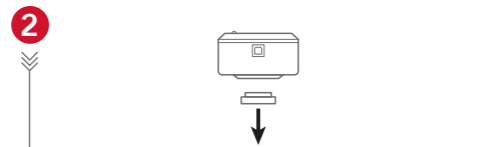
ES Accesorios incluidos en el paquete:
DE Im Paket enthaltenes Zubehör:
FR Accessoires inclus dans l'emballage :
IT Accessori inclusi nella confezione:

1. ES Moticam
DE Moticam
FR Moticam
IT Moticam
2. ES Aro C
DE C-Ring
FR Anneau C
IT Anello a C
3. ES Objetivo enfocable (12 mm)
DE Fokussierbare Linse (12 mm)
FR Objectif focalisable (12 mm)
IT Lente foceggiabile (12 mm)
4. ES Adaptadores de ocular (2)
DE Okularadapter (2)
FR Adaptateurs pour oculaires (2)
IT Adattatori per oculare (2)
5. ES Cable USB
DE USB-Kabel
FR Câble USB
IT Cavo USB
6. ES Macro-tubo
DE Makrotubus
FR Macro-tube
IT Tubo macro
7. ES Preparación de calibración
DE Kalibriermastab
FR Lame de calibrage
IT Vetrino di calibrazione
8. ES Software (enlace de descarga)
DE Software (Link zum Herunterladen)
FR Logiciel (lien de téléchargement)
IT Software (link per il download)

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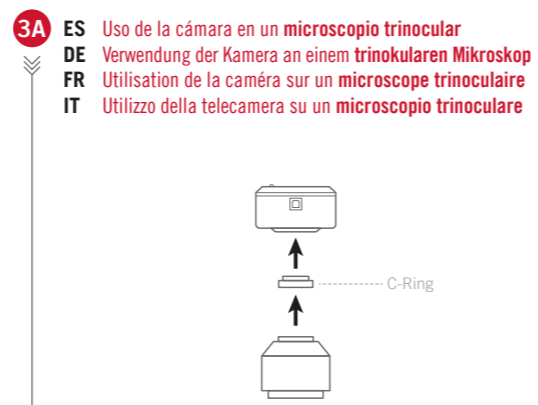
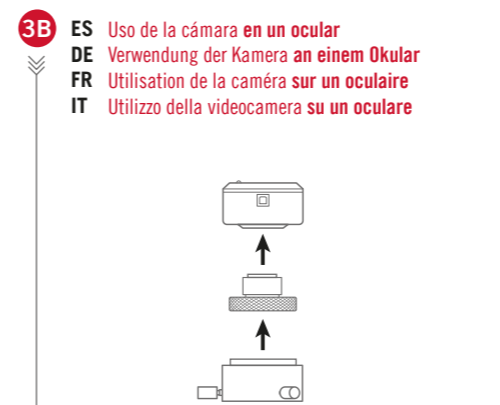
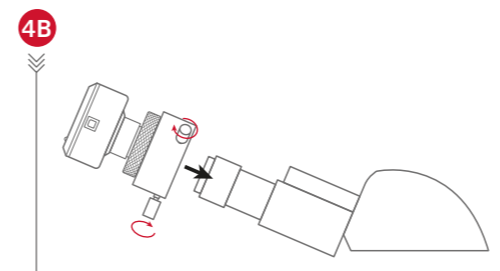
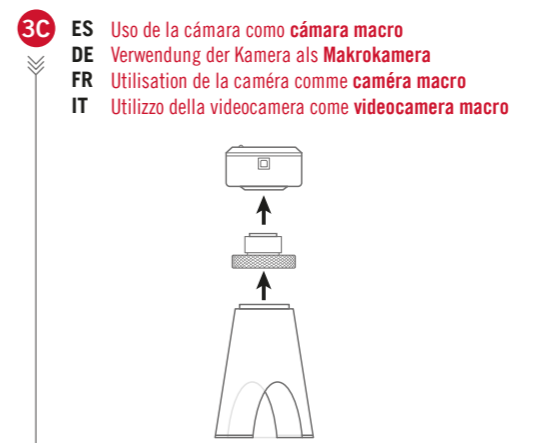
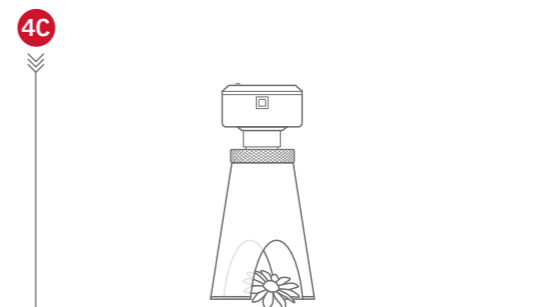
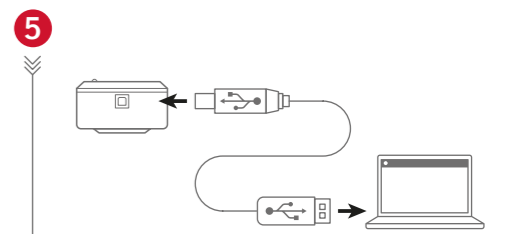
MotiConnect **FIRST STEPS**

MotiConnect **CALIBRATION & MEASUREMENTS**

North America

support@motic-america.com
1-877-901-4141

Europe

digital.support@moticeurope.com
+34 93 756 62 86

Global

info_digital@motic.com

EN This camera is a delicate instrument: please handle it with care and, when not in use, store it in a safe place with the dust cap attached. Avoid contact with water and exposing the device to extreme temperatures (below 0 or above 50 degrees Celsius). To allow us to offer you the highest level of service, please retain your receipt as proof of purchase and register your camera online at www.motic.com today. Our customer service team is at your disposal to assist you with any queries.

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Motic Asia

Motic Headquarters (Hong Kong)

Unit 2002, L20, Tower Two, Enterprise Square Five,
38 Wang Chiu Road, Kowloon Bay, Kowloon
Tel: 852-2837 0888 | Fax: 852-2882 2792

Motic Xiamen (China)

4A-6, N°810 Fangshan Nan Road, XiangAn,
Xiamen, P.R. China.
Tel: 0592 559 9681 | Fax: 0592 559 9683

Design Change: The manufacturer reserves the right to make changes in instrument design in accordance with scientific and mechanical progress, without notice and without obligation.

*CCIS® is a trademark of Motic Incorporation Ltd. Motic Incorporation Limited Copyright © 2002-2023. All Rights Reserved. ISO 9001:2000, ISO 14001:2004 and ISO 13485 Certified.

Created and designed in Barcelona (Spain) by Motic Europe | May 2023