



# PIG® Hazardous and Chemotherapy Drug Spill Kit – Low Volume

## Includes:

- 2 - PIG® High-Visibility Absorbent Mat Pads (absorbs up to 832 mL)
- 1 - Protective Apron with Sleeves (Polyethylene Coated Polypropylene)
- 1 - Pair of XL Boot Covers (Polyethylene Coated Polypropylene)
- 2 - Pairs of Size Large Nitrile Gloves (ASTM D6978)
- 1 - Chemotherapy Waste Disposal Bag (Yellow 20-gallon)

## Instructions:

### Step 1 — Wear appropriate PPE

Put on appropriate personal protective equipment (PPE) prior to entering the spill area.

### Step 2 — Absorb the liquid

Working from the outside of the spill inward, place mats on the spill until the liquid is completely covered and absorbed.

### Step 3 — Dispose of used absorbents

Place all used absorbents or other contaminated materials in the disposal bags and tie the bag closed. Dispose of the bag and all used materials in compliance with local, state and federal regulations.

## Kit antiderrames PIG para sustancias peligrosas y fármacos de quimioterapia - Volumen bajo

## Incluye:

- 2 tapetes absorbentes PIG® de alta visibilidad (absorben hasta 832 ml)
- 1 mandil protector con mangas (polipropileno recubierto con polietileno)
- 1 par de cubiertas extragrandes para botas (polipropileno recubierto con polietileno)
- 2 pares de guantes de nitrilo talla grande (ASTM D6978)
- 1 bolsa para desechos de quimioterapia (color amarillo, 20 galones)

## Instrucciones:

### Paso 1 — Utilice el EPP adecuado




Utilice el equipo de protección personal (EPP) adecuado antes de ingresar al área del derrame.

### Paso 2 — Absorba el líquido

Coloque los tapetes sobre el derrame, desde la parte exterior y trabajando hacia adentro, hasta que el líquido quede cubierto y se absorba completamente.

### Paso 3 — Deseche los absorbentes usados

Coloque todos los absorbentes utilizados, así como otros materiales contaminados, en las bolsas de desechos y amárrelas para cerrarlas. Deseche la bolsa y todos los materiales utilizados de modo que cumpla con las reglas locales, estatales y federales.

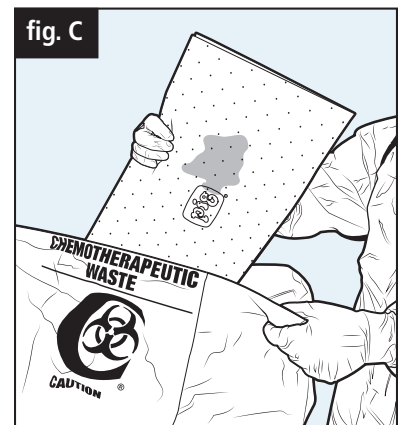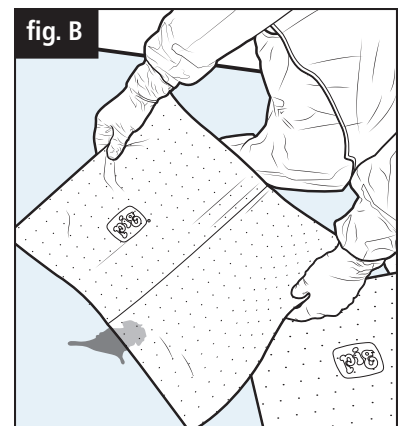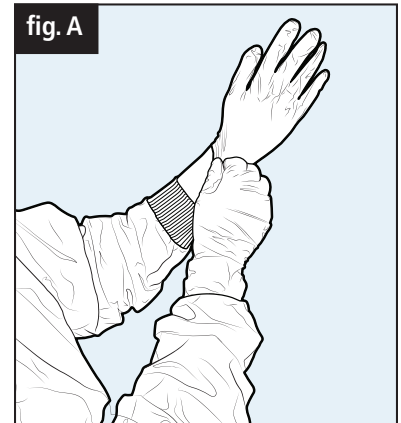

**New Pig**

One Pork Avenue  
Tipton, PA 16684-0304

[newpig.com](http://newpig.com)

North America: **1-800-468-4647**

Europe: **+31 (0)76 596 92 50**

China: **+86-21-400 921 5178**

PIG, PIG logo are registered trademarks in USA and other countries. See [tm.newpig.com](http://tm.newpig.com)

UK: **0800 919 900**

Outside North America: **+1-814-684-0101**

166071 / 358751